



GeneQuery™ HLA-A PCR+Sanger SBT Typing Kit (SBT-A)

Catalog #GK891
40 samples

Product Description

The human leukocyte antigen (HLA) complex is a gene complex located on the short arm of human chromosome 6. The HLA complex is the human version of the major histocompatibility complex (MHC) which helps the immune system distinguish host cells from invaders. It is the most polymorphic locus in the human genome, and has been implicated in many pathologies including autoimmunity disorders and cancer.

GeneQuery™ HLA-A PCR+Sanger SBT Typing Kit (SBT-A) utilizes the Sequence-Based Typing (SBT) method, which is the gold standard for HLA typing. This method involves PCR amplification of HLA genes, followed by Sanger sequencing of the amplicons. The acquired sequences are aligned to an HLA allele database using the SOAPTyping software (BGI Genomics) to obtain the 4-digit genotypes of the samples. The two haplotypes from both parents can be typed simultaneously using this kit and the SOAPTyping software.

SBT-A kit contains a ScienCell proprietary HLA-A amplification primer set (A_ampF primer mix and A_ampR primer mix; Cat #GK891a) that can amplify a region in the human genome that covers exons 2-4 of HLA-A gene, and 6 sequencing primer mixes (A_2F, A_2R, A_3F, A_3R, A_4F, and A_4R; Cat #s GK891b-g) that can be used in the following Sanger sequencing of the PCR product. Please refer to Figure 2 for the binding location of the primers on the human genome. Each GeneQuery™ HLA typing kit contains a human Genomic DNA Reference (GDR) sample with known HLA alleles (Cat #GK891r). The GDR can be used to monitor the quality of the reagents and the instruments. The HLA-A alleles in GK891r are HLA-A*02:01 and HLA-A*25:01.


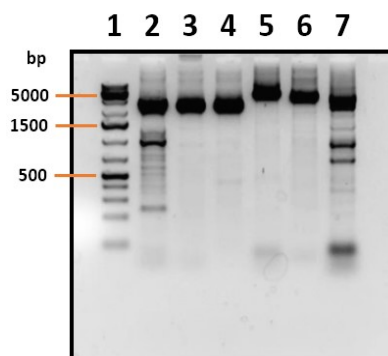


Figure 1. Gel electrophoresis separation of PCR amplification products with human genomic DNA as the template using ScienCell GeneQuery™ HLA PCR+Sanger SBT Typing Kits.

Lane 1, DNA ladder;
Lane 2, PCR product of HLA-A amplification PCR primer set (Cat #GK891a, size: ~3,100 bp);
Lane 3, PCR product of HLA-B amplification PCR primer set (Cat #GK892a, size: ~2,800 bp);
Lane 4, PCR product of HLA-C amplification PCR primer set (Cat #GK893a, size: ~2,900 bp);
Lane 5, PCR product of HLA-DPB1 amplification PCR primer set (Cat #GK894a, size: ~4,900 bp);
Lane 6, PCR product of HLA-DQB1 amplification PCR primer set (Cat #GK895a, size: ~4,300 bp);
Lane 7, PCR product of HLA-DRB1 amplification PCR primer set (Cat #GK896a, size: ~3,600 bp);

Figure 2. Primer binding locations on the human genome.

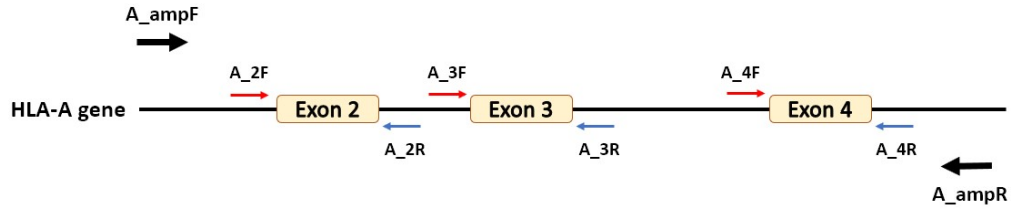

Kit Components

Cat #	Component	Quantity	Storage
GK891a	HLA-A amplification primer set (A_ampF and A_ampR)	400 µL	-20°C
GK891b	A_2F sequencing primer mix, 10 µM	200 µL	-20°C
GK891c	A_2R sequencing primer mix, 10 µM	200 µL	-20°C
GK891d	A_3F sequencing primer mix, 10 µM	200 µL	-20°C
GK891e	A_3R sequencing primer mix, 10 µM	200 µL	-20°C
GK891f	A_4F sequencing primer mix, 10 µM	200 µL	-20°C
GK891g	A_4R sequencing primer mix, 10 µM	200 µL	-20°C
GK891h	Nuclease-free H ₂ O	4 mL	4°C
GK891r	Human genomic DNA reference (GDR), ~50 ng/µL (HLA-A*02:01 and HLA-A*25:01)	10 µL	-20°C
MB6038-1	2X EmeraldNStart HiFi Marathon PCR Master Mix, 1 mL	2 vials	-20°C

Additional Materials Required (Materials Not Included in Kit)

Component	Recommended
genomic DNA template	Customers' samples
PCR purification kit	SpeedDNA PCR Purification Kit (ScienCell, Cat #MB6948)
Sanger sequencing kit	
PCR plates or tubes	

Quality Control

The primers are validated by PCR, followed by gel electrophoresis and Sanger sequencing using GK891r as the template.

Product Use

SBT-A is for research use only. It is not approved for human or animal use, or for application in clinical or *in vitro* diagnostic procedures.

Shipping and Storage

SBT-A is shipped on dry ice. Upon receipt, store the primers and EmeraldNStart HiFi Marathon PCR Master Mix at -20°C in a manual defrost freezer, and nuclease-free H₂O at 4°C. Once the EmeraldNStart HiFi Marathon PCR Master Mix is thawed, store it at 4°C and do not refreeze. It is stable at 4°C for up to 6 months if stored properly.

Procedures

Important: *Only use nuclease-free reagents in PCR amplification.*

1. Thaw the HLA-A amplification primer set (Cat #GK891a). Aliquot as needed. Store at -20°C in a manual defrost freezer. Avoid repeated freeze-and-thaw cycles.
2. For each genomic DNA sample, prepare one 100 µl PCR reaction solution as shown in Table 2. Aliquoting to multiple PCR tubes for the PCR reaction to keep the volume in each tube low is recommended.

Table 2.

Genomic DNA template	80 - 120 ng
HLA-A amplification primer set (Cat #GK891a)	10 µl
2X EmeraldNStart HiFi Marathon PCR Master Mix	50 µl
Nuclease-free H ₂ O (Cat #GK891h)	variable
Total volume	100 µl

3. Seal the PCR reaction tubes. Centrifuge the tubes at 1,500x g for 15 seconds.
4. Refer to Table 3 for PCR program setup.

Table 3.

Step	Temperature	Time	Number of cycles
Initial denaturation	95°C	10 min	1
Denaturation	95°C	30 sec	38
Annealing	65°C	30 sec	
Extension	72°C	4 min	
Final extension	72°C	10 min	

5. Refer to the Instructions for Use of the PCR purification kit to be used for the purification of the PCR products. Purified PCR products can be used for the Sanger sequencing.
6. Refer to the Instructions for Use of the Sanger sequencing kit to be used for the sequencing of the PCR products. Six sequencing reactions should be prepared with each of the 6 sequencing primers included in the kit.
7. Analyze the sequencing results with the SOAPTyping software (BGI Genomics) to obtain the 4-digit genotypes of the samples. The software can be downloaded at <https://github.com/BGI-flexlab/SOAPTyping/releases> (for Windows, Linux, and Mac) or <https://www.sciencellonline.com/products-services/genetics-genomics/soaptyping> (for Windows only). Please refer to <https://github.com/BGI-flexlab/SOAPTyping> for Instruction for Use of the SOAPTyping software. Appendix 1 provides a quick start guide for the SOAPTyping software. If allele-specific sequencing primers (ASSPs) are required to resolve the typing ambiguity, please refer to Appendix 2 for ASSPs designing guide.

Appendix 1. Quick start guide for the SOAPTyping software

Disclaimer: SOAPTyping Software is a free and open-source software developed by BGI Genomics. ScienCell is NOT responsible for technical support of the software.

1. SOAPTyping


The screenshot displays the SOAPTyping software interface with several key components highlighted by red boxes and labels:

- Toolbar:** Located at the top left, containing various icons for file operations and analysis.
- Base Navigator:** Located below the toolbar, showing a genomic map with Exon2, Exon3, and Exon4, and a sequence alignment view.
- Sequence Display:** A detailed view of the DNA sequence with a yellow highlight on a specific region.
- Chromatogram Display:** A chromatogram showing the signal intensity for each base (A, C, G, T) corresponding to the sequence above.
- Sample/file List:** A list of sample files on the right side, including 1462_A_2F_A05.ab1, 1462_A_2R_D05.a_2R, 1462_A_3F_A01.ab1, 1462_A_3R_F05.ab1, 1462_A_4F_C05.ab1, and 1462_A_4R_G05.ab1.
- Allele Match List:** A table at the bottom right showing allele matches for different samples, with columns for Allele1, Allele2, Info, and GSSP.
- Nucleotide Codes:** A bar at the bottom left showing the sequence R-AGY-TCK-GTM-ACS-GCW-ATB-GTCD-GATH-ACTV-GCA.
- Base Information:** A bar at the bottom right showing Exon:2 Codon:-114, 2 Pos:269 Code:AAW QV:30.


- Verify that the names of the .ab1 files to be analyzed are in the format of “Sample name”_“HLA gene-of-interest”_“Exon targeted and primer direction”_“Additional info”.ab1. “HLA gene-of-interest” is “A” when using this kit for typing HLA-A gene. “Additional info” is a user-defined annotation to the file. The SOAPTyping software uses the underscore symbol “_” to separate the sections. It is required to have 4 sections in the file name and no additional underscore symbol in the file names. For example,

- 1462_A_2F_A05.ab1
- 1462_A_2R_D05.ab1
- 1462_A_3F_B05.ab1
- 1462_A_3R_F05.ab1
- 1462_A_4F_C05.ab1
- 1462_A_4R_G05.ab1

3. Run the SOAPTyping software, click “Open File”, then click “Open” button to select files for analysis.

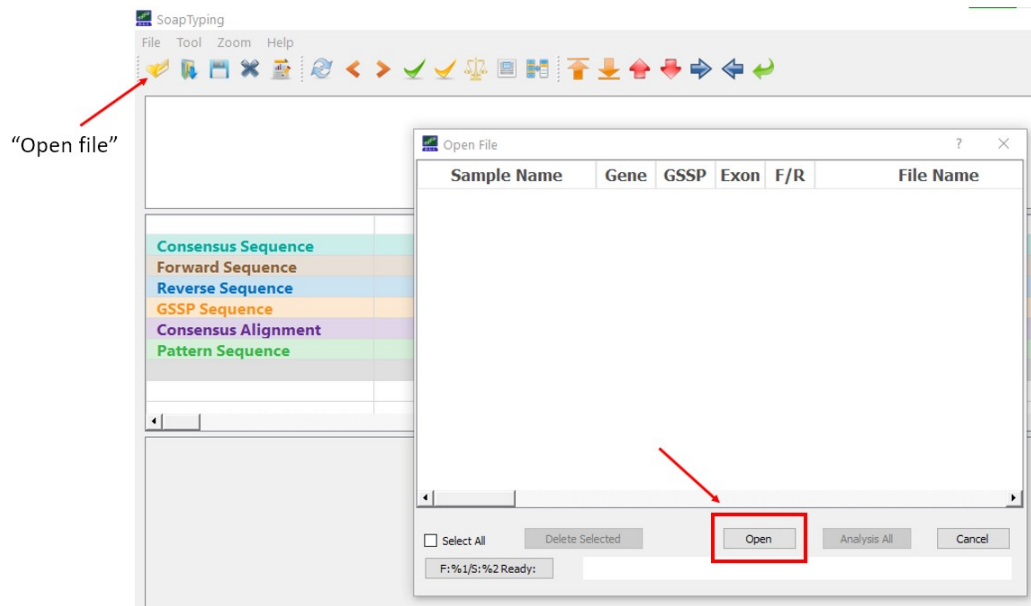

4. Select and open the files, then click “Analyze all” button to analyze the data. The “Gene”, “Exon”, and “F/R” sections should be filled automatically. If not, please check if the file names follow the format described in Step 2 of Appendix 1.

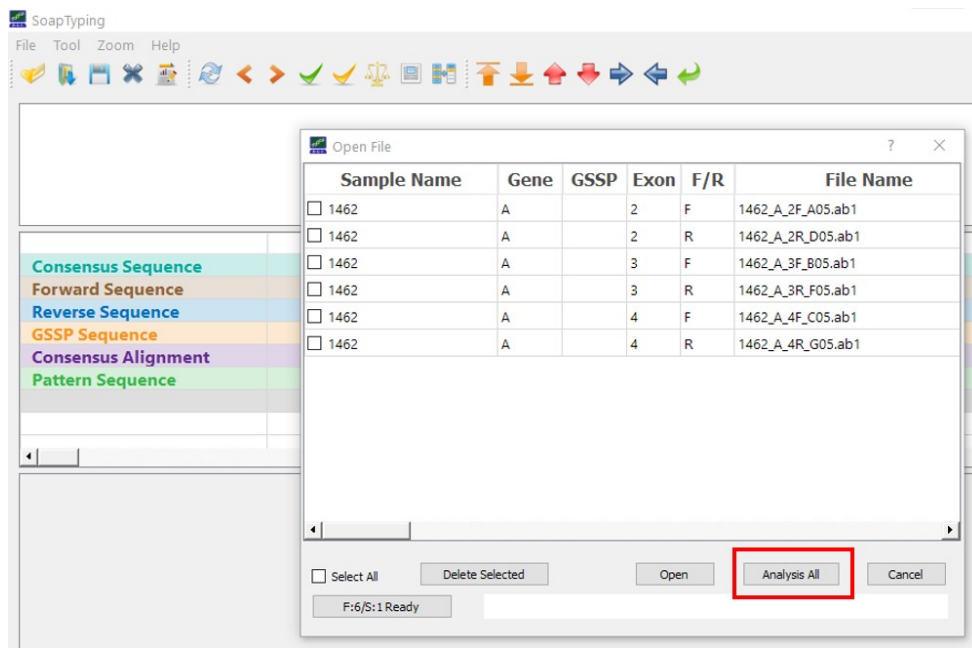


5. The results with ambiguities and mismatches to resolve will be listed.




In the upper half of the Base Navigator panel, the ambiguities and mismatches are indicated by color-coded vertical lines. Blue lines indicate different but compatible forward and reverse sequences (e.g., W and A); Green lines indicate different and



The peak of T was NOT accounted for, therefore A should be W here

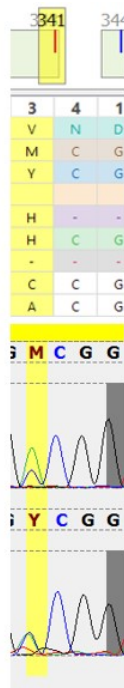

The peak of T should not be picked, therefore Y should be M here

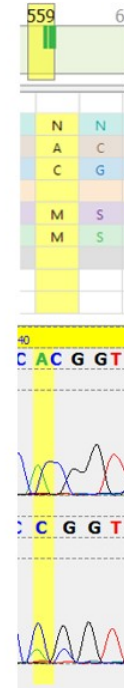

This position should be an M. In forward sequencing, A and C were separated. In reverse sequencing the peak of A was NOT accounted for, causing a mismatch in forward and reverse sequences.

incompatible forward and reverse sequences (e.g. S and T); Red lines indicate differences between consensus sequence from database and pattern sequence of the sample; Black lines indicate manually changed and now compatible nucleotides.

The red lines in the lower half of the Base Navigator panel indicate mismatches between the pattern sequence and the sequence of the selected allele pair from the Allele Match List panel. These ambiguities and mismatches are usually caused by the interpretation of the Sanger sequencing results. All the ambiguities and mismatches in green and red must be manually corrected to get the final types approved.



